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

1. Remove the microscope cover.
2. Turn on the PC (①). Log in automatically (User name: Leica-User).  
※If user name is "LASX-Admin", change it to "Leica-User" from the lower left corner of the screen and log in. User name: **Leica-User**, Password : **none**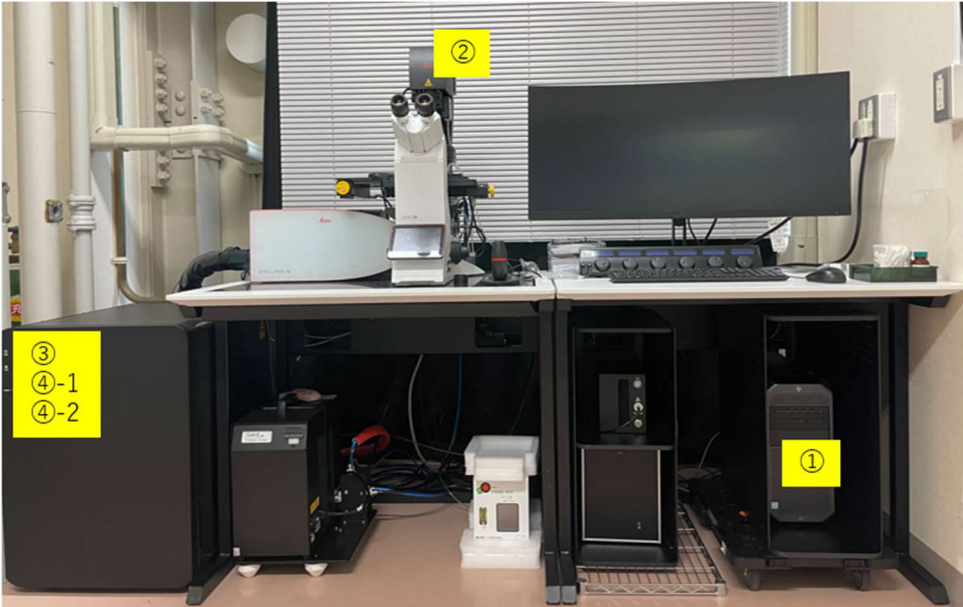

3. Tilt the condenser unit (②) of the microscope backward.

**【caution】**

Check that the gas supply tube extending to the rear of the stage is as shown in the figure below.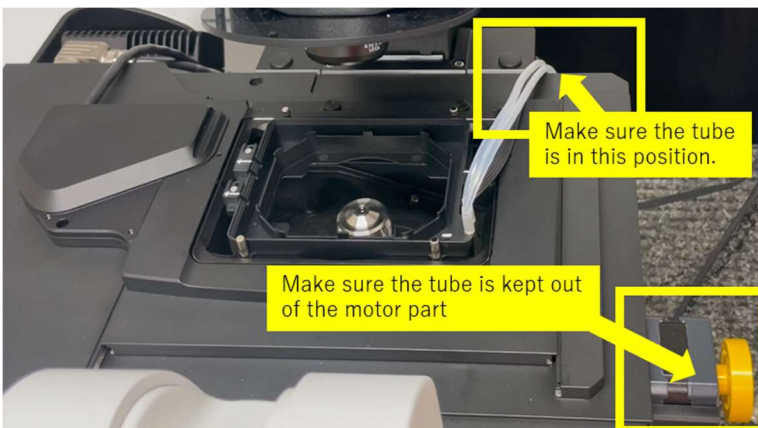

4. Turn on main switch (③).


**【caution】** The stage initialization begins.

**Make sure the tube is kept out of the motor part!**

After initialization completed, put the condenser unit (②) back into place.

5. Turn on the laser (④-1).

6. Turn the key switch (④-2) 90° clockwise.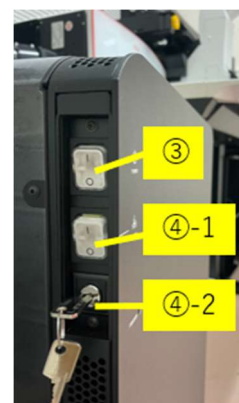

7. Double-click the LAS X software shortcut icon on the desktop to launch the software.

Select options as below and click OK.

Configuration: **machine.xlhw**, Microscope: **DMI8**, Load settings at startup: **OFF**


1. Select 10x objective lens.

※Make sure 40x and 60x cleaned properly if you used!

2. Turn off the laser on the software.

Click “Configuration” tab, and select “Laser Config”. Turns off all lasers.

**【Caution】** Do not change the WLL output %. (85% is the default at our facility)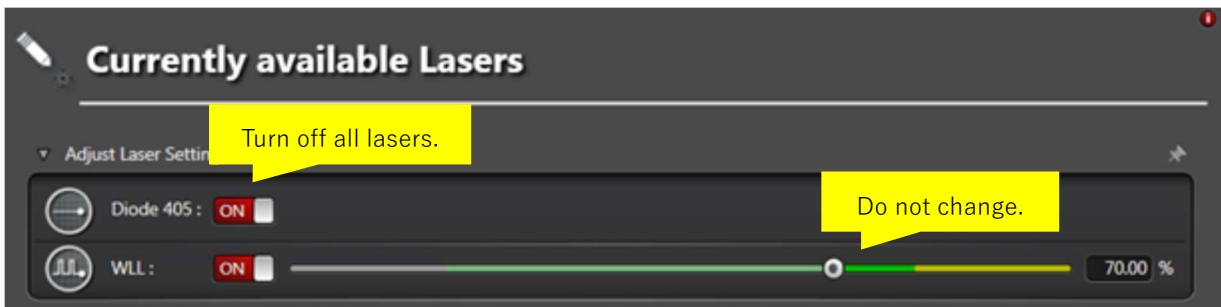


3. Close LAS X software.

4. Turn off the PC.

5. Turn the key switch (④-2) 90° counterclockwise.

6. Turn off the laser (④-1).

7. Turn off main switch (③).

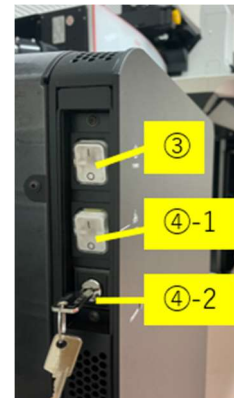### Use only LAS X software for data analysis

1. Turn on the PC (①). Log in automatically (User name: Leica-User).

2. Double-click the LAS X software shortcut icon on the PC to launch the software.  
Select options as below and click OK.

Configuration: **Simulator.xlhw**, Microscope: **DM Manual-6**, Load settings at startup: **OFF**

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