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XGEN CUSTOM HYB PANEL, XGEN MRD PANEL

INTEGRATED DNA TECHNOLOGIES

For research use only. Not for use in diagnostic procedures. Unless otherwise agreed to in writing, IDT does not intend these products to be used in clinical applications and does not warrant their fitness or suitability for any clinical diagnostic use. Purchaser is solely responsible for all decisions regarding the use of these products and any associated regulatory or legal obligations. The information in this presentation is confidential and proprietary. No part of the presentation may be used, reproduced or distributed without the express written permission of Integrated DNA Technologies, Inc.

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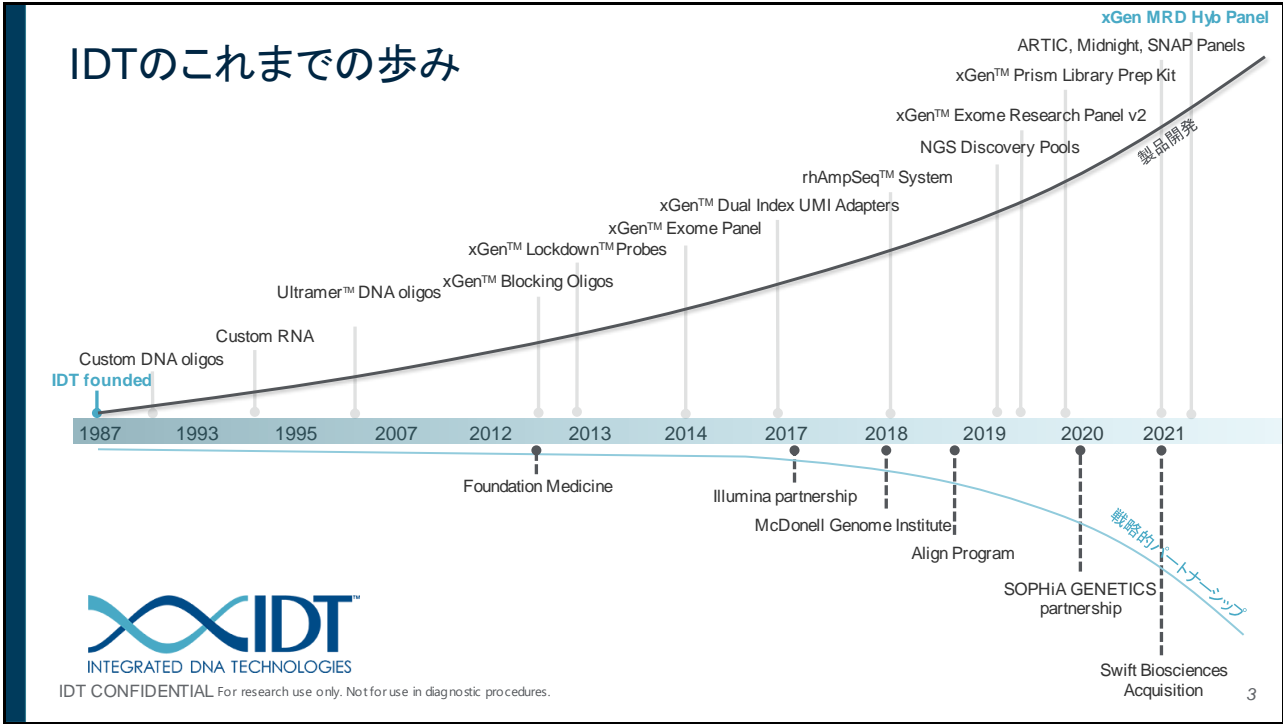
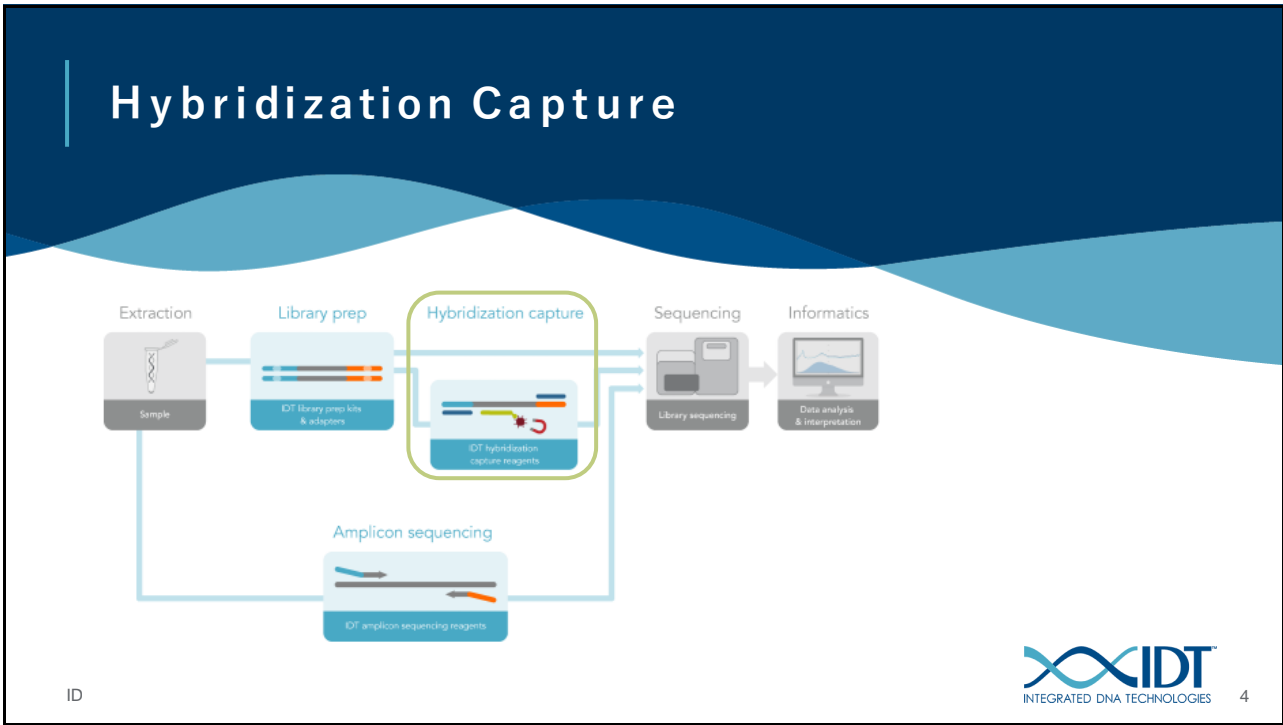
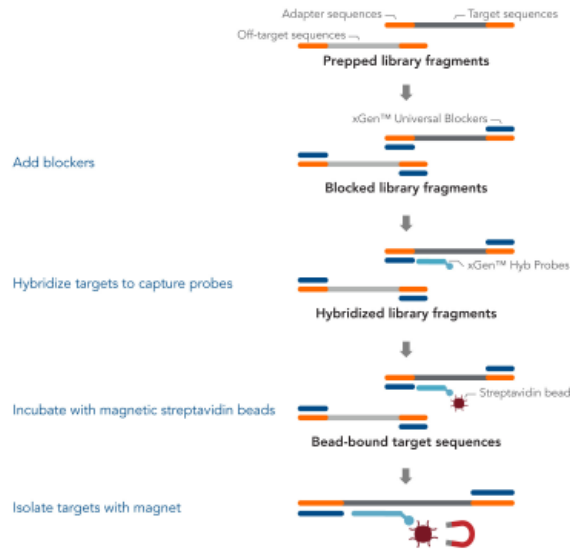
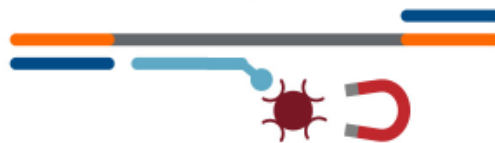
6

xGen Hybridization Capture Core Reagents

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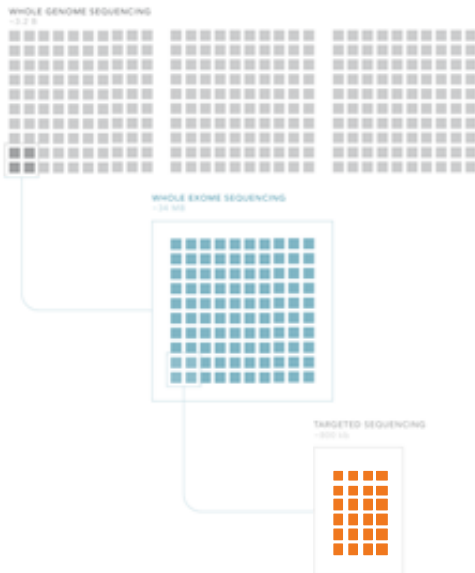
Flow Cytometry

Next Generation Sequencing (NGS)

For research use only. Not for use in diagnostic procedures.

11

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with our xGen WGS Solution

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with our xGen WES or other Targeted Solutions

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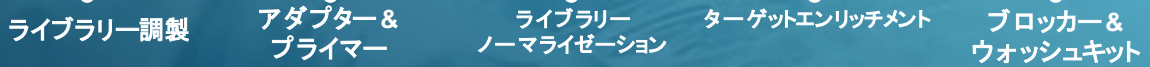
with our xGen MRD Solution

For research use only. Not for use in diagnostic procedures.

IDT CONFIDENTIAL

12


xGen - NGS made for MRD research


IDT CONFIDENTIAL For research use only. Not for use in diagnostic procedures.

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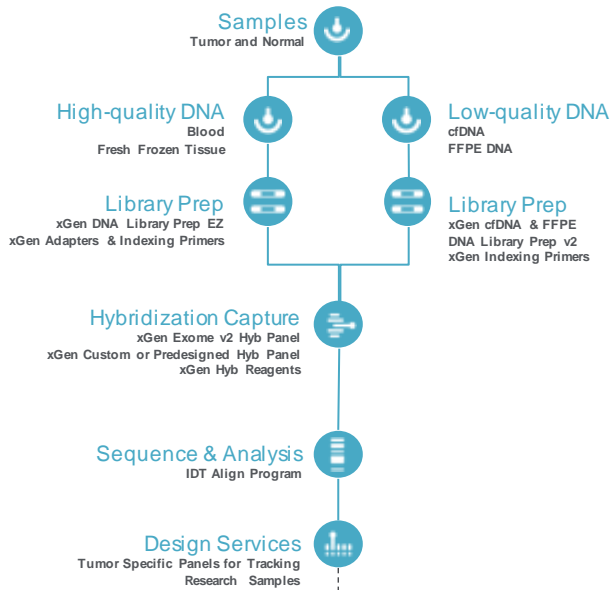
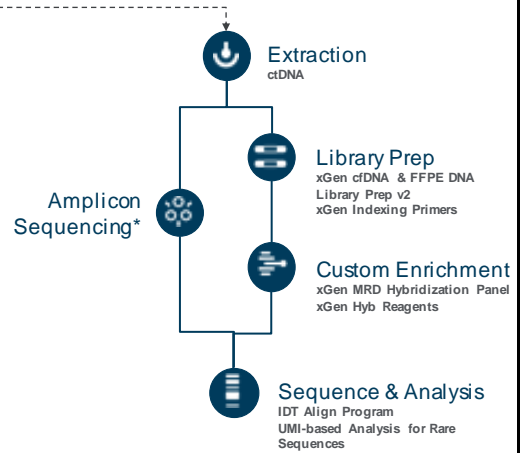
For research use only. Not for use in diagnostic procedures.

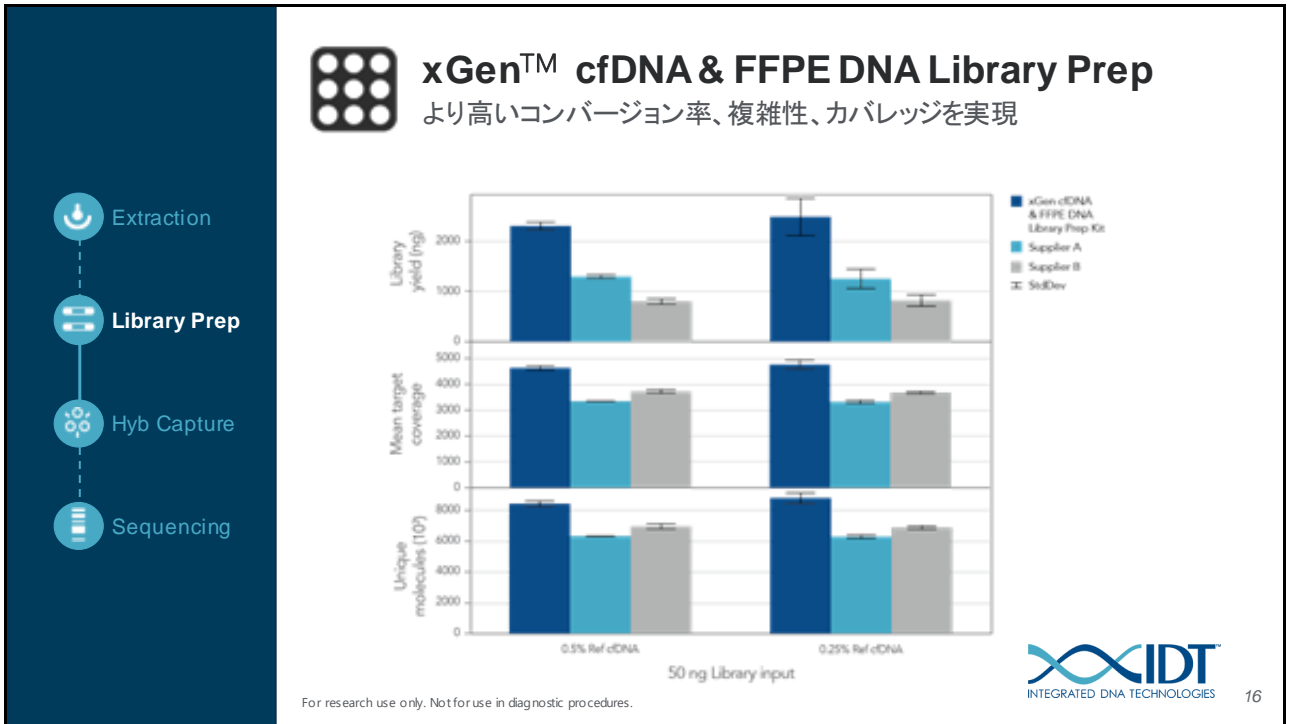*consult your IDT sales representative to learn more about IDT's amplicon sequencing solutions 14

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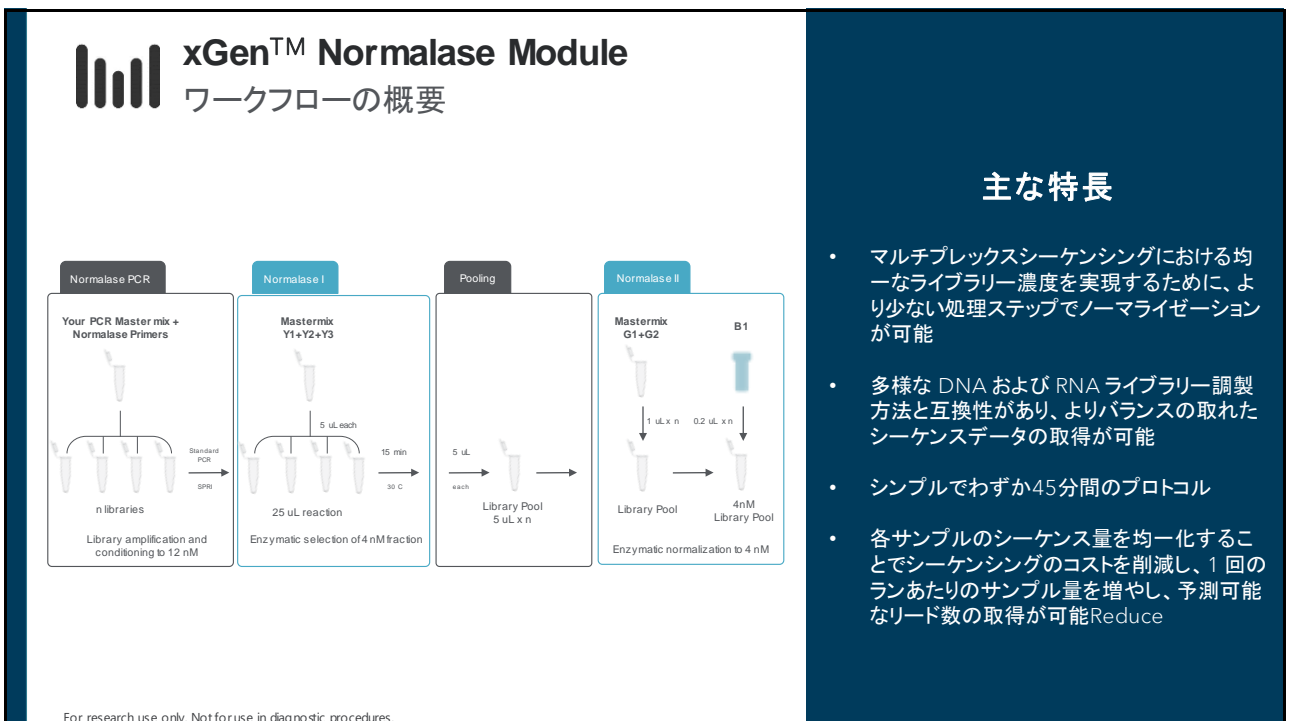Feature	xGen Custom Hyb Panel—Accel	xGen Custom Hyb Panel	xGen Custom Hyb Panel—Production (No functional testing)	xGen Custom Hyb Panel—Production
Custom design service	✓	✓	✓	✓
Quality synthesis	✓	✓	✓	✓
NGS Functional Test Report		✓		✓
Quality documentation			✓	✓
Stock custom panel	✓	✓	✓	✓
Reaction sizes	16, 96, custom	16, 96, custom	16, 96, custom	16, 96, custom
Turnaround time	~2週間	通常~4週間	Project dependent	Project dependent

23

| Custom Hyb Panel定価価格表 – 16 rxn

Max # of Probes	16 rxn			
	xGen Custom Hyb Panel - Accel	xGen Custom Hyb Panel	xGen Custom Hyb Panel - Production No Functional Test	xGen Custom Hyb Panel - Production
500	¥290,000	¥395,000	¥527,200	¥595,000
1,000	¥350,000	¥440,000	¥580,000	¥810,000
5,000	¥585,000	¥990,000	¥1,398,100	¥1,445,000
7,500	¥655,000	¥1,110,000	¥1,530,000	¥1,810,000
10,000	¥700,000	¥1,145,000	¥1,810,000	¥2,210,000
25,000	¥1,750,000	¥1,985,000	¥4,520,000	¥4,880,000
50,000	¥3,500,000	¥3,735,000	¥9,045,000	¥9,175,000

24

| Custom Hyb Panel定価価格表 – 96 rxn

Max # of Probes	96 rxn			
	xGen Custom Hyb Panel - Accel	xGen Custom Hyb Panel	xGen Custom Hyb Panel - Production No Functional Test	xGen Custom Hyb Panel - Production
500	¥875,000	¥1,005,000	¥1,355,100	¥1,620,000
1,000	¥1,165,000	¥1,285,000	¥1,985,000	¥2,135,000
5,000	¥1,750,000	¥2,160,000	¥2,805,000	¥3,200,000
7,500	¥2,190,000	¥2,550,000	¥3,280,000	¥3,920,000
10,000	¥2,390,000	¥2,735,000	¥3,350,000	¥4,070,000
25,000	¥2,915,000	¥3,210,000	¥4,665,000	¥5,155,000
50,000	¥3,790,000	¥3,995,000	¥9,335,000	¥9,825,000

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0 ITEMS \$0.00 USD

[Order by stock part number](#)

PRODUCTS & SERVICES ▾ APPLICATIONS & SOLUTIONS ▾ SUPPORT & EDUCATION ▾ TOOLS ▾ COMPANY ▾

xGen™ Hyb Panel Design Tool

An updated solution offering enhanced design functionality to generate your custom hybridization capture panels for targeted sequencing on Illumina platforms.

- Easy design for even coverage across targets
- Capable of generating targeted sequencing panel designs for multiple genomes with custom tiling options
- Fast, expert consultation available from our technical support team to address custom specifications outside the scope of the publicly available tool

Sign In

Keep me signed in. [Details](#) [Forgot Password](#)

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xGen Hyb Panel Design Tool

Welcome to the IDT xGen Hyb Panel design tool. If you would like a consultation to address a design outside the scope of this tool, submit a request to the xGen custom design service here.

Create new design Dashboard

Input format: BED target coordinates

Species: Homo sapiens (Human) NCBI GRCh37.p13 (hg19) - 105.20201022

Run probe QC: Yes No

Panel name: Name your design...

Probe length: 120

Probe tiling density: 1x

Paste/Type input: Upload file

Submit your BED coordinates here. Both .txt and .bed file extensions are acceptable.
Example: BED File

Choose File NO FILE CHOSEN

DESIGN
CLEAR AND RESET
Need design assistance?

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Design Tool Dashboard

My xGen Designs

Panel name	Source	Date & Time	Probe count	Info	ACTIONS
Elephant Gene target 1	IDT Design	6/25/2021, 3:27:23 PM	2805	In process	ACTIONS
Human Gene Target 1	Design Tool	6/25/2021, 2:46:45 PM	2942	In process	ACTIONS

Show/Hide Inactive Designs

Need design assistance? Contact ngsdesign@idtdna.com or have our Design Team assist with a Custom Design.

Design date / time	Species	Target base count	Probe Count	Percent covered	Output Files
6/25/2021, 2:46:45 PM	Homo sapiens (Human) NCBI GRCh38.p13 (hg38)	45668	2942	99.2	Download

Quote suggestions

xGen Custom Hyb Panel

xGen Custom Hyb Panel - Accel

xGen Custom Hyb Panel - Production

xGen Custom Hyb Panel - Production (No functional testing)

Need more options? Select "Custom"

16 rxn

96 rxn

Custom rxn

Comments & Customizations

Specify desired panel customizations here for our Design Team.

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xGen™ MRD Hybridization Panel Ordering Tool

Configure

Selected product: xGen™ MRD-Hyb Panel, 24 rxn

Minimum probes per panel: 50

Maximum probes per panel: 2,000

Probe length restriction: 60 - 120

Maximum panels: 50

Supported file format: Excel spreadsheet (.xls, .xlsx) **Download Template**

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	A	B
1	Panel name	Probe Sequence
2	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
3	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
4	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
5	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
6	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
7	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
8	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
9	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
10	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
11	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
12	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
13	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
14	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
15	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
16	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
17	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
18	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
19	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
20	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
21	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
22	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
23	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG
24	lipid_panel_x1	TGCTCCAGGCGGGGCGGAAATTTTCTGAGGGATTTTGTACTGCGCGGGCGGCGGGGCGG

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- xGen Broad-range RNA Library Prep Kit
- xGen RNA Library Prep Kit

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- xGen Methyl-Seq DNA Library Prep Kit

38

THANK YOU

39